


Guard Column Holder for 1 cm Guard Columns - Manual (P/N 424267)

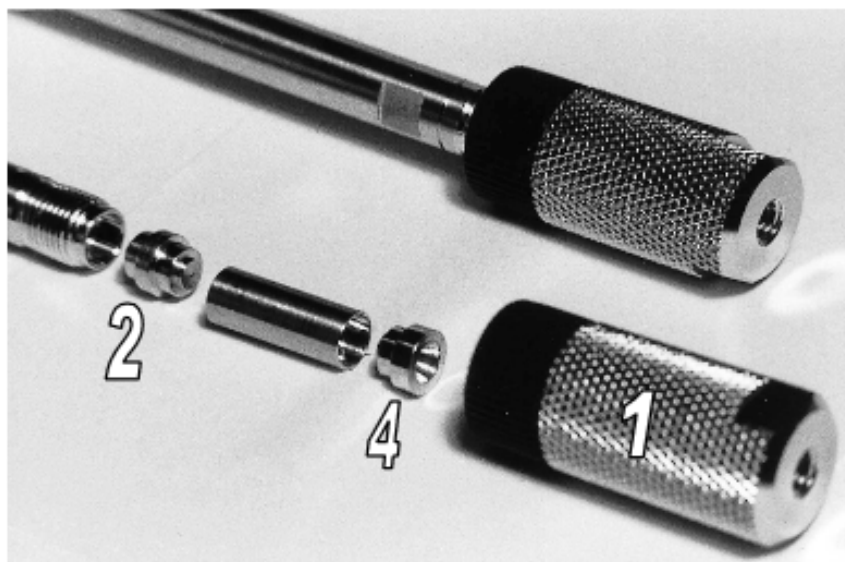

The guard column holder set contains: guard column holder (1)
connector (2)

For usage of the guard column holder the column head (3) of the main column and the fitting adapter (4) which is placed at the column entrance are being removed (if needed with a gripper)

Within the column the micronic sieves with overlying PTFE seal are now visible. The stainless steel connector (2) needs to be placed right upon that seal.

Remove the protecting caps from the 1 cm guard column and put the guard column upon the connector (2). Place the afore removed fitting adapter (4) on the other side of the guard column.

Put the guard column holder (1) over the guard column incl. fitting adapter and tighten by hand (if needed with a wrench).

The solvent capillary can now be screwed into the inner thread (thread: UNF 10/32 for 1/16" screw ports) of the guard column holder.